



Instruction for Use

AcuQ 1X dsDNA HS Assay Kit

AG12549

Version.V1E1

**Research Use Only
Not For Diagnosis Procedures**

1. Description

This product is a fluorescent quantification kit designed for rapid, sensitive, and accurate detection of double-stranded DNA (dsDNA) when used with Qubit fluorometers. The kit includes a premixed working solution (containing fluorescent dye) and dsDNA standards. In a 200 µl reaction system, it can accurately quantify dsDNA samples within a concentration range of 10 pg/µl to 100 ng/µl (with sample volumes of 1–20 µl, corresponding to a total dsDNA amount of 0.2–100 ng). The assay demonstrates strong tolerance to common contaminants such as proteins, salts, and detergents within a defined range. Compared to single-stranded DNA (ssDNA), RNA, proteins, and free nucleotides, this kit exhibits high selectivity for dsDNA.

The kit features a ready-to-use premixed formulation for simple and convenient operation. To perform the assay, simply mix the target dsDNA sample with the working solution, incubate at room temperature for 2 minutes, and measure using a Qubit fluorometer.

2. Kit Information

Kit Name	Cat. No	Specification
<i>AcuQ</i> 1X dsDNA HS Assay Kit	AG 12549	100 Rxns

3. Transportation and Storage

Storage	Store at 4°C
Transportation	Transport at 4°C / Ice Bag

4. Kit Components

Kit Components	Volume
1X dsDNA HS Working Solution	50 ml
dsDNA Standard 1	1 ml
dsDNA Standard 2	1 ml

5. Protocol

1. Pre-experiment Preparation

- ① Before use, equilibrate all kit components to room temperature.
- ② Prepare an adequate number of Qubit assay tubes and label them appropriately. Avoid labeling on the sidewalls of the tubes to prevent interference with fluorescence signal detection.


2. Sample Preparation

- ① Standard preparation: Take two 0.5 mL Qubit assay tubes, add 190 µl of 1X dsDNA HS Working Solution to each, followed by 10 µl of dsDNA Standard 1 and dsDNA Standard 2, respectively. Vortex gently for 2-3 sec to mix, avoiding bubble formation, and briefly centrifuge.
- ② Test sample preparation: Add 180-199 µl of 1X dsDNA HS Working Solution to a 0.5 mL Qubit assay tube, then introduce 1-20 µl of the test sample to achieve a final volume of 200 µl per tube. Vortex gently for 2-3 sec to mix, minimize bubbles, and centrifuge briefly.

3. Detection

- ① Incubate all Qubit assay tubes at room temperature for 2 min, protected from light.
- ② Follow the Qubit fluorometer operating instructions and select the 1X dsDNA High Sensitivity assay program for measurement.

Experimental Workflow

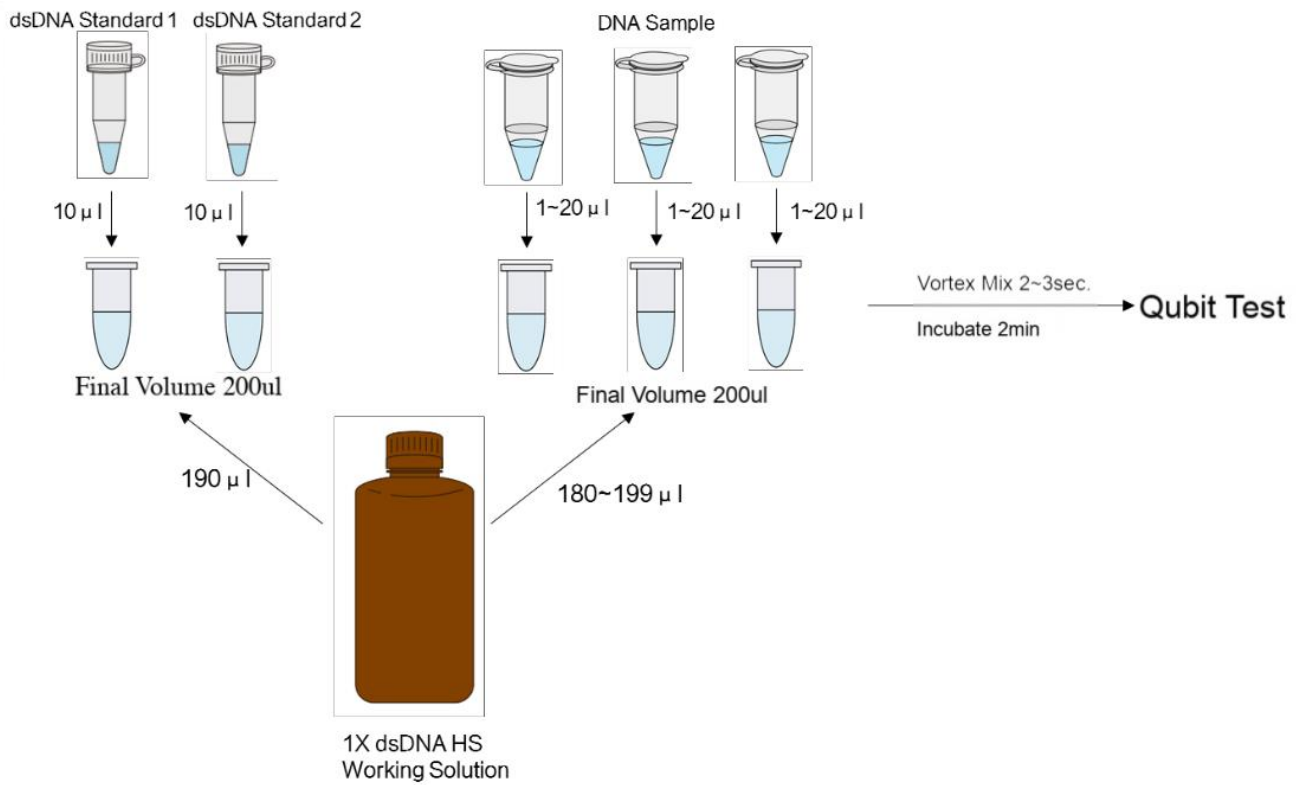

Accurate Biotechnology (Hunan) Co., Ltd

No.1 JinYang Road, WangCheng District, Changsha City, Hunan Province, China

service@agbio.com.cn

+86 400 767 6022

en.agbio.com.cn

Research Use Only